# Software for Spectroradiometer

## CS-S30

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English

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## Software License Agreement

The terms of the license agreement of this software are provided in the [Software License Agreement] dialog box displayed on-screen during the installation process. This software can be installed only if you agree to all the terms of the agreement.

### Safety Precautions

Before using the software, please read this manual as well as the instruction manuals of your instrument and personal computer (PC) carefully to ensure safe and correct use.

Every effort has been made to ensure the accuracy of the content of this software. However, should you have any questions or comments, please contact the nearest KONICA MINOLTA authorized service facility.

## Notes on Use

- This software is designed to be used with the Windows 10, Windows 11, macOS Catalina, macOS Big Sur, macOS Monterey, macOS Ventura operating system. Note that no operating system is included with this software.
- One of these operating systems must be installed on the PC before this software can be installed.

## About the Instruction Manual

- The instruction manual will be saved to the PC as a PDF file during installation of the software. Foreign-language versions of the instruction manual PDF file will also be saved.

## Formal designations of application software used in this manual

(Designation in this manual)	(Formal designation)
Windows, Windows 10	Microsoft® Windows® 10 Pro Operating System
Windows, Windows 11	Microsoft® Windows® 11 Pro Operating System
macOS, macOS Catalina	Apple® macOS® Catalina
macOS, macOS Big Sur	Apple® macOS® Big Sur
macOS, macOS Monterey	Apple® macOS® Monterey
macOS, macOS Ventura	Apple® macOS® Ventura

## Trademarks

- Microsoft, Windows, Windows 10, and Windows 11 are registered trademarks of Microsoft Corporation in the U.S.A. and other countries.
- Apple, macOS Catalina, macOS Big Sur, macOS Monterey, macOS Ventura are registered trademarks of Apple Inc. in the U.S.A. and other countries.
- Intel and Core are registered trademarks of Intel Corporation in the U.S.A. and other countries.
- Other company names and product names mentioned in this manual are the registered trademarks or trademarks of their respective companies.

# 1. System Environment

## Operating requirements

<b>OS</b>	Windows 10 Pro 64-bit(×64) Windows 11 Pro macOS Catalina macOS Big Sur macOS Monterey macOS Ventura (Rosetta 2 is required when using a Mac with an Apple Silicon chip installed.)
	<ul style="list-style-type: none"> <li>• The required PC system configuration is the recommended configuration for the operating system above or the specifications below (whichever is more advanced).</li> </ul>
<b>CPU</b>	At least as advanced as the Intel® Core™ i Series At least as advanced as the Apple Silicon M1 chip
<b>Memory</b>	4 GB or more
<b>Hard disk</b>	At least 100 MB of free space. At least 50 MB of the hard disk's free space needs to be on the system drive (the drive where the operating system is installed).
<b>Display resolution</b>	Must support at least 1,280 × 768 pixel, 16 bit color display
<b>Other</b>	USB port supporting at least USB 2.0 is needed for instrument connection USB port supporting at least USB 2.0 is needed to attach the protection key (when connecting and using the CS-2000 or CS-2000A)

# 2. Installing/Uninstalling the Software

The software can be downloaded from the following website.

<https://www.konicaminolta.com/instruments/download/software/display/index.html>

## 2.1 Using Windows

Install the CS-S30 software.

- Be sure to log in to the PC as a user with Administrator privileges to install this software.
- A User Account Control dialog will be displayed several times during installation of the software and the drivers. Click [Continue] for each dialog.

### 2.1.1 Installing the Software

1. Start Windows.
2. Download the installer (cs-s30\_verxxx\_win\_multi.zip) from the URL above. (“xxx” indicates the version number.)
3. Unzip the downloaded ZIP file. Verify the destination folder for the extracted files.
4. Double-click the extracted setup file (setup.exe). The installation program will start.
  - The installation program will examine the PC being used. It may take several minutes for this process to be completed and for the setup dialog to be displayed.
5. A dialog box for selecting the language to use during setup will be displayed.
  - Select the language to use during installation. To select a different language, click on the currently selected language and select the desired language from the displayed pull-down menu.
6. The installation program may notify you of the need to install Microsoft components. If this occurs, the installation of the necessary components will begin automatically.
  - Agreement to the licenses for the necessary components being installed will be required.
  - The PC may need to be restarted. In such cases, restart the PC and proceed from Step 1 above.
  - Installation may take a long time (10 minutes or more).
7. Read through the software license agreement and select “Accept” and “Next” to agree and continue with the installation.
8. Verify the installation destination folder. To use the selected folder for installation, click [Next].
  - To change the destination folder, click [Browse...], specify the desired folder, and then click [OK].
9. Select whether to create desktop icons using the appropriate checkbox. Click [Next].
10. Verify the settings and click [Install] to begin installation.
  - Installation may take several minutes. Do not perform any other operations on the PC during installation.
11. Finalize installation. Click [Finish].

### 2.1.2 Uninstalling the Software

1. Right-click the Start button on the screen.
2. Select [Apps and Features] from the menu.
3. Select [CS-S30 Verxxx] and click the [Uninstall] button. (“xxx” indicates the version number.)
4. A dialog box confirming uninstallation of the software will be displayed. Click [Uninstall] to proceed.
5. A dialog box confirming removal of the software will be displayed. Follow the on-screen instructions to uninstall the software.

## 2.2 Using macOS

### 2.2.1 Installing the Software

1. Download the installer (cs-s30\_verxxx\_mac\_multi.dmg) from the URL above. (“xxx” indicates the version number.)
2. Double-click the downloaded DMG file. The CS-S30 drive will be displayed.
3. Double-click “setup.pkg.” The installation program will start.
4. Click [Continue].
5. Read through the license agreement and click [Continue]. A confirmation dialog box will be displayed. Click [Agree] to agree to the conditions and continue with the installation.
6. Set the destination for installing the software. Click [Change Install Location...] to change the destination as necessary.
7. Click [Install]. Copying of the files will begin. This may take several minutes. You will be asked to enter the password for the PC to permit installation.
8. Finalize installation. Click [Close].

When starting this software for the first time, a screen that guides you to install Rosetta may appear.

In that case, click [Install] to install Rosetta. During the process, you will be asked to enter a password to authorize the installation.

After installing Rosetta, start this software again.

### 2.2.2 Uninstalling the Software

Drag the CS-S30 app from the [Applications] folder to the Trash.

## 3. Connecting

Use a USB cable to connect the PC to the spectroradiometer. Refer to the appropriate instruction manual for more information on hardware-based connections. Refer to the instruction manual for this software for more information on configuring the software connection settings.

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